# SUPAPORE NP



### Pleated Nylon Membrane Cartridge

SupaPore NP Microfiltration cartridges use a pleated high surface area Nylon 66 membrane that has been designed for use in a wide range of liquid processing applications. The naturally hydrophilic Nylon 66 membrane contains no wetting agents and has very low levels of extractables making this cartridge an ideal choice for systems requiring a reliable and consistent high-quality filtrate.

#### **Product Features**

SupaPore NP cartridges feature:

- A naturally hydrophilic N66 membrane
- Exceptionally high flow rates with low clean pressure losses
- Wide chemical compatibility

All cartridges are thermally bonded and adhesive free. Each unit is preflushed with ultra pure water and integrity tested before final assembly.

- Non Pyrogenic
- Materials meet US FDA Title 21 and USP Class VI requirements
- Meet current USP requirements for extractables, particle release and endotoxins
- Fast rinse-up time to 18 Megohm-cm\*1
- Available in a wide range of absolute rated pore sizes.

Suitable for most sanitisation regimes including steaming, autoclaving, hot water flush and most common sanitising agents.

(Note: Steaming, autoclaving and hot water sanitisation must use the Reinforced Polypropylene End Cap option)

Amazon Filters manufactures a comprehensive range of filter housings for SupaPore NP cartridges including the industrial 50 and 60 Series and the 70 Series for hygienic applications. Please consult our Sales Office for further details on the full range.



### Features and Benefits

- High surface area giving high flow rates and low initial pressure losses
- Wide chemical compatibility
- Meets FDA requirements
- Thermal bonding process eliminates adhesives and ensures minimal extractables
- Identification mark on every cartridge
- Comprehensive Product Validation Guide available
- All individual packing bags are anti static

### Industries and Applications

Pharmaceuticals Fine Chemicals Ophthalmics, Oral medications, Bioburden reduction

Fine Chemicals
Electronics

Aqueous and Organic bulk feeds, Solvents, Bubble jet inks High purity water, Polishing loop return lines, Point of use

Cosmetics

Alcohol, Mouthwashes, Make up water



### SupaPore NP Technical Data

#### **Dimensions**

Outside Diameter: 68.5mm
Typical Surface Area: 0.65m² (10")

#### Sterilisation and Sanitisation\*3

 $\begin{array}{lll} \mbox{Steam or Autoclave:} & 121^{\circ}\mbox{C for 15 mins (40 cycles)} \\ \mbox{Hot Water:} & 90^{\circ}\mbox{C for 30 mins (0.2 bar $\Delta p$ max)} \\ \mbox{$^{*3}$ Applies to single open end cartridges only. For all steaming and hot water applications, the Reinforced Polypropylene End Cap option must be used} \end{array}$ 

#### **Maximum Operating Conditions**

Temperature: 80°C

Recommended Maximum Differential Pressure: Forward Flow: 4.0 Bar @ 20°C Reverse Flow: 3.5 Bar @ 20°C


Recommended change-out differential pressure: 2.5 Bar

Product Validation Guide available on request. All **SupaPore NP** cartridges are manufactured under strict control with batch number identification, giving full traceability on all components.

#### **Standard Materials of Construction**

Filter Media: Nylon (N)
Media support: Polypropylene
End Caps: Polypropylene
Cage/Core: Polypropylene

#### Flow Rates (10" Element)

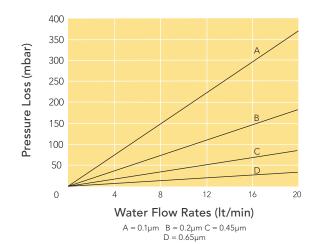

## Amazon Filters Ltd.

Albany Park Estate, Camberley, Surrey, GU16 7PG, ENGLAND

Tel: +44 (0) 1276 670 600 Email: sales@amazonfilters.co.uk Web: www.amazonfilters.com